


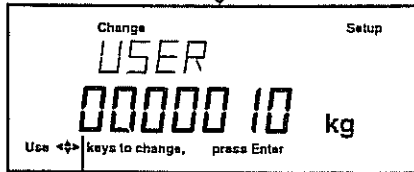
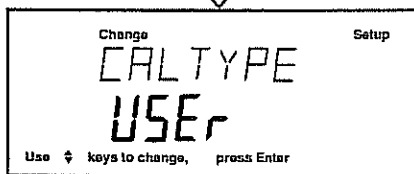


### 4.1.2 User Calibration

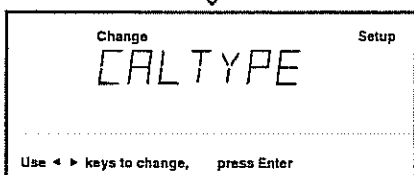

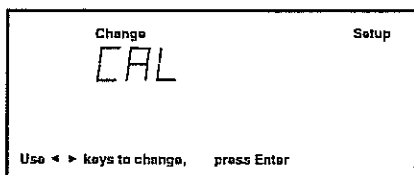

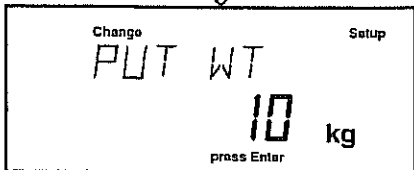

User calibration is used when it is desired to calibrate the Indicator using a mass of known value from 10% to 100% of the capacity of the scale base. To use this calibration feature, proceed as follows:



(FLASHING)



### Procedure

- Press the **Setup** button, SCALE is displayed.
- Press  button UNTIL CAL is displayed.
- Press **Enter** button, CAL TYPE is displayed.
- Press **Enter** button, SPAN is displayed.
- Press  button to select user calibration. CAL TYPE USEr is displayed.
- Press **Enter** button; the display indicates the last calibration mass value which was entered with the first digit flashing. (Sample illustrates 10kg).
- Press   and or   and enter the desired mass value when using Model CD-31. The numerical keypad can be used for Model CD-33. This number must be at least 10% of the full span value.
- Press **Enter** button, CLR PAN is displayed.
- Press **Enter** button, WORKING is momentarily displayed, then PUT WT 10 kg is displayed.
- Place specified calibration mass on pan.
- Press **Enter** button, WORKING is displayed. After a few seconds CAL SET is displayed, the display then returns to WEIGH mode and indicates mass on pan.
- Remove calibration mass from the pan.
- User calibration is completed.





Ohaus Corporation  
29 Hanover Road  
Florham Park, NJ  
07932-0900

# INSTRUCTION MANUAL

## *Models CD-31 and CD-33 Indicators*





↓  
indicator  
CD-33 35AS

mark correction program

1000

MA 2000 kg

1000 10 07000

Attn: Repair Dept

Program 1000  
unit

Large Pan

Small Pan

Wt = Kg

Wt = ~~2%~~ Kg

Zero = 2%

Zero = 2%

Cal unit = Kg

Cal unit = Kg

Grad = 0.001

Grad = 0.001

Full SC = 7 digits 50 Kg

Full SC = 7 digits 10 Kg

Lock = off

Lock = off

Ohaus Corporation  
29 Hanover Road,  
Florham Park, NJ 07932, USA  
Tel: (973) 377-9000,  
Fax: (973) 593-0359

With offices in England, Germany, France, Spain, Italy, Poland, Mexico, Japan, Korea and China.



15331700A